

Dharmacon™

RNAi, Gene Expression & Gene Editing

DharmaFECT™ Duo Transfection Reagent – for siRNA and Plasmid Co-Transfection

Introduction:

Dharmacon DharmaFECT Duo is a lipid based reagent specially formulated for co-transfection of plasmid and siRNA. Under optimized conditions, efficient delivery of both plasmid and siRNA can be achieved with minimal cell toxicity. As is generally observed with plasmid transfection, cell viability is reduced compared to siRNA transfection alone. When evaluating delivery conditions, a threshold of 70% cell viability and 80% gene silencing is set for successful, non-toxic co-transfection of plasmid and siRNA.

Product Description:

- Provided as a liquid at a concentration of 1 mg/mL.
- Confirmed to be free of microbial contaminants and ribonucleases.

Product	Size	Cat. #
DharmaFECT Duo	0.2 mL	T-2010-01
	0.75 mL	T-2010-02
	1.5 mL	T-2010-03
	1.5 mL × 5 tubes	T-2010-04

Shipping and Storage:

- DharmaFECT Duo co-transfection reagent is shipped on ice packs. If the ice melts during transit there is no risk to the product. Stability testing demonstrates no loss in product functionality after prolonged storage at warm temperatures (see the Stability Testing Product Bulletin for additional details).
- Upon receipt, store at 4 °C. **DO NOT FREEZE.** DharmaFECT transfection reagents are labeled with a date of manufacture and are good for 24 months following that date when stored as recommended.

Related Products:

- We recommend including a positive and negative control, such as Dharmacon RNAi Control Reagents, in every RNAi experiment. For more information, click [here](#).

Accompanying Documents:

- siRNA/Plasmid Co-Transfection Protocol

Publication Reference Guide:

When referencing the use of DharmaFECT Duo Co-Transfection Reagents, please include the following information: DharmaFECT Duo Transfection Reagent, GE Healthcare Dharmacon, Inc., Lafayette, CO.

GE, imagination at work and GE monogram are trademarks of General Electric Company. Dharmacon is a trademark GE Healthcare companies. All other trademarks are the property of General Electric Company or one of its subsidiaries. ©2014 General Electric Company—All rights reserved. Version published June 2014. GE Healthcare UK Limited, Amersham Place, Little Chalfont, Buckinghamshire, HP7 9NA, UK



GE Healthcare

Orders can be placed at:
gelifesciences.com/dharmacon

Customer Support: cs.dharmacon@ge.com
Technical Support: ts.dharmacon@ge.com or
1.800.235.9880; 303.604.9499 if you have any questions.